


# Sampling in an Aseptic Process

## Risk Mitigation & Regulatory Compliance



Somasundaram G.  
Senior Technical Consultant, Asia Pacific  
Process Solution, Merck KGaA  
October 12, 2020

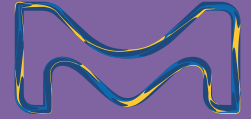

The life science business  
of Merck KGaA, Darmstadt, Germany  
operates as MilliporeSigma in the  
U.S. and Canada

**Millipore®**


Preparation, Separation,  
Filtration & Monitoring Products

# Agenda

- 1 Introduction to Sampling – Where, What & Challenges
- 2 Regulatory recommendations and corresponding needs – Key Drivers
- 3 Complexity and Risks of traditional sampling methods
- 4 Key considerations for Closed & Disposable sampling options

# Introduction to Sampling Where, What & Challenges

## Aseptic processing – A path full of pitfalls...

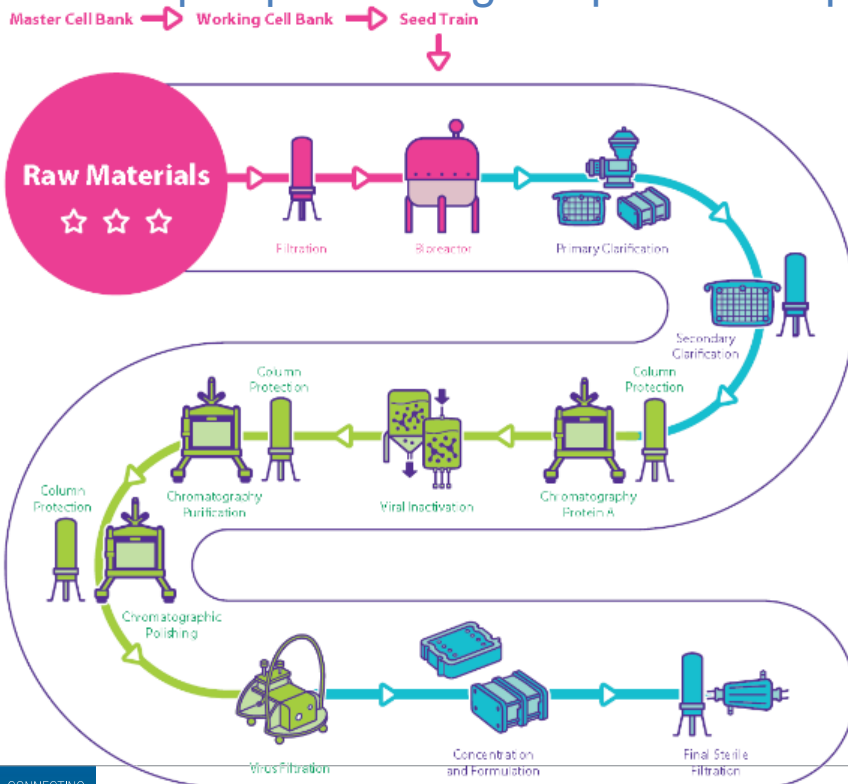


Biopharma manufacturing operations require substantial precautions to prevent microbial ingress:

- 1 Highly trained personnel
- 2 A risk-based approach to contamination control and a dedicated strategy
- 3 Resulting in significant costs:
  - Infrastructure
  - Day-to-day operations

# Introduction to sampling

## What are The Requirements and Challenges?

- Representative Sample :
  - Composition
  - Non-operator dependent
  - Repeatable
- Protect the Process (Contamination-free)
- Protect the Person (Contained)
- Protect the Product (Contained)
- Easy of Use
- Save Time and Cost Efficient

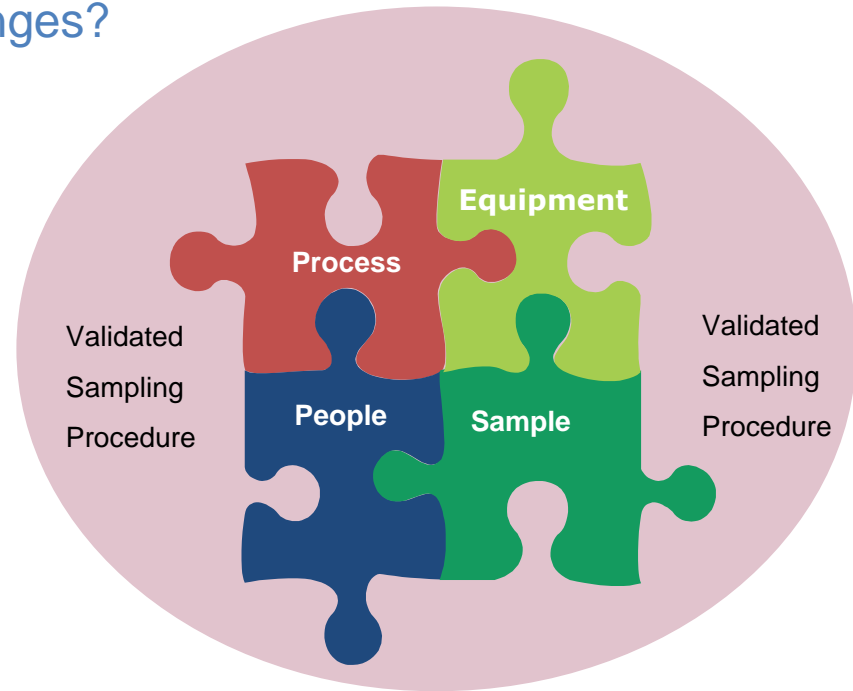

In today's biopharmaceutical market, sampling is critical during every step in the manufacturing process. An imprecise or false positive result can lead to a quarantine as well as the need to repeat the analysis.

# Sampling in Bioprocesses

	Bioreactor Seed train	Buffer/Media Preparation	Bioreactor Production	Purification	Sterile Filtration
pH	✓	✓	✓	✓	✓
Conductivity/ Osmol	✓	✓	✓	✓	✓
Cells	✓		✓		
Metabolites	✓		✓		
Protein analysis			✓	✓	✓
Bioburden & Archiving	✓	✓	✓	✓	✓
Endotoxin	✓	✓	✓	✓	✓
Other (virus, by products)	✓	✓	✓	✓	✓


# Sampling plan & options

## Final Filling


Where to Sample



Sampling point

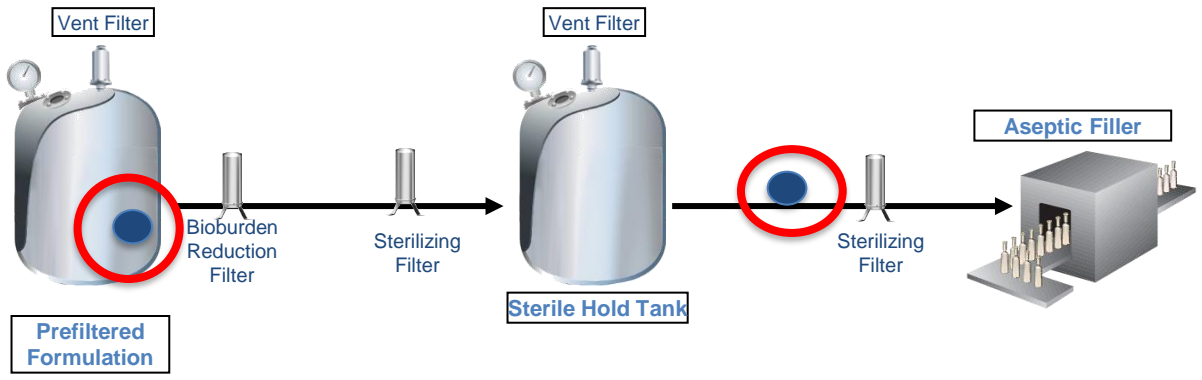

Single Sterilizing Filter

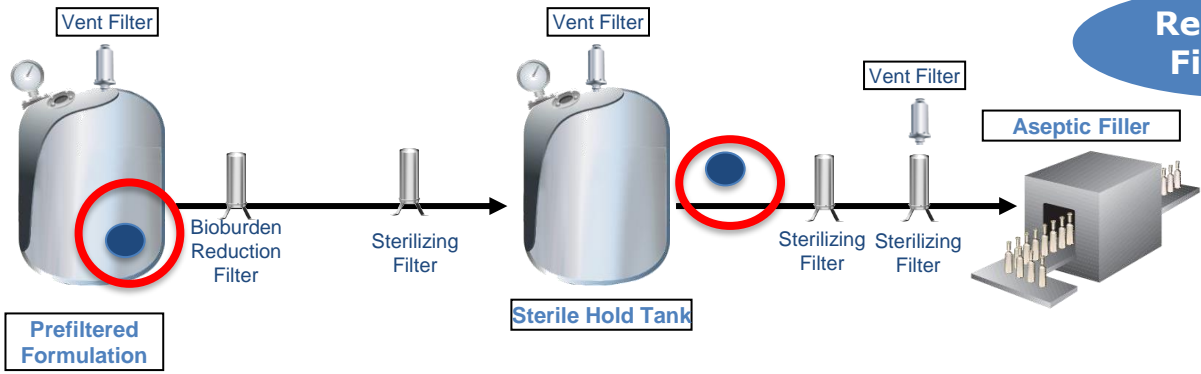


Redundant Filtration





# Regulatory recommendations and corresponding needs – Key Drivers

# Drivers for Aseptic Process Sampling

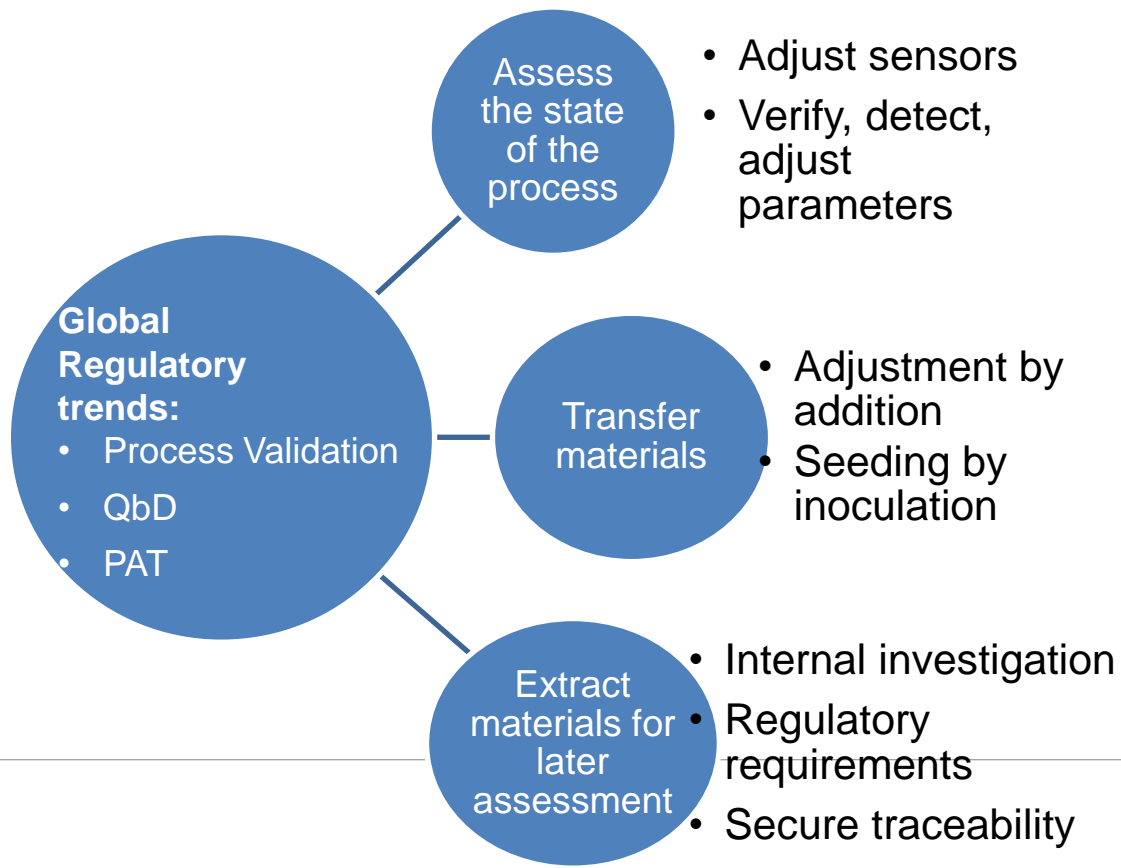

# Relevant Regulatory Agencies & Industry Associations

Regulatory Agencies / Industry Associations	Relevant Documentation
Food and Drug Administration	cGMP guidance
World Health Organization	<b>Annex 2</b> – Good manufacturing guidance for API <b>Annex 4</b> - Good Manufacturing Practices for pharmaceutical products: main principles
European Medicines Agency	GMP Annex 1 - Manufacture of Sterile Medicinal Products
International Council for Harmonization of Technical Requirements for Pharmaceuticals for Human Use	ICH Q7 – GMP guidance for API
Parenteral Drug Association	Report #69 - Bioburden and Biofilm Management in Pharmaceutical Operations
The Pharmaceutical Inspection Convention and Pharmaceutical Inspection Co-operation Scheme	Guide to good manufacturing practice for medicinal products – Part I

# Extracts of Relevant Requirements

Recommendations	Source
Contamination control & Monitoring before bioburden reduction	<ul style="list-style-type: none"> <li>• <b>PICS/S – FDA cGMP</b> – Part 1 §5.19 –f               <ul style="list-style-type: none"> <li>• Use of closed system recommended from phase 1</li> </ul> </li> <li>• <b>WHO Annex 4</b> <ul style="list-style-type: none"> <li>• “The use of disposable sampling materials has distinct advantages”</li> </ul> </li> <li>• <b>WHO Annex 2 - ICH Q7A</b> - GMP guidance for API</li> <li>• <b>EU GMP Annex 1</b></li> <li>• <b>EudraLex</b> – Vol 4 – Part II – 2009</li> </ul>
Operator bias elimination	<ul style="list-style-type: none"> <li>• <b>WHO Annex 4</b></li> <li>• <b>FDA</b></li> <li>• <b>European Pharmacopeia</b> - Guidelines for Sampling of Pharmaceutical Products and Related Materials</li> </ul>
Representative sample	
Health & safety focus	
Retained samples	<ul style="list-style-type: none"> <li>• <b>FDA cGMP</b> Guidance for the industry investigational drugs section F. Laboratory Controls / 1. Testing</li> <li>• 2 years after expiration date / completion of trial and twice the quantity necessary to perform all tests</li> </ul>

# Sampling Process & Precautions

## WHO Annex 4: Preparation for sampling

- “All sampling tools and implements should be made of inert materials and kept scrupulously clean. [...]”
- “The cleaning procedure used for all sampling tools and implements should be documented and recorded.”
- “The use of disposable sampling material has distinct advantages”

## WHO Annex 4: Sampling Operations and Precaution

“There should be a written procedure describing the sampling operation. [...]. It should ensure that representative samples are taken in sufficient quantity for testing in accordance with specifications.”

## WHO Annex 4: Storage and Retention

“The container used to store a sample should not interact with the sampled material nor allow contamination. [...]. As a general rule the container should be sealed and preferably tamper-evident.”



**World Health  
Organization**



# Sampling of Pharmaceutical Products and Related Materials



- FDA cGMP for phase 1 drugs  
“recommends the use of closed system to minimize the risk of contamination”
- Guidance for the industry investigational drugs section  
F. Laboratory Controls / 1. *Testing*  
“We recommend that the sample consist of a quantity adequate to perform additional testing or investigation if required at a later date [...]. We recommend that you appropriately store and retain the samples for at least two years [...].”
- Q7A GMP guidance for manufacturing API – section C. In-process Sampling and Controls (8.3)  
“In-process sampling should be conducted using procedures designed to prevent contamination [...]. [...] ensure the integrity of samples after collection.”

# Annex 1 EU GMP (Draft)

## Regulatory requirements

**8.94 Bioburden samples should be taken from the bulk product and immediately prior to the final sterile filtration. Systems for taking samples should be designed so as not to introduce contamination.**

**9.7 Sampling methods should not pose a risk of contamination to the manufacturing operations.**


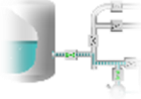






**7.3 Non-essential processes such as product inspection and in process testing should be conducted with the crimping option there is a validated and easy disconnection in place to further treat the sample outside the clean areas wherever possible.**

# Complexity & Risk of Traditional Sampling Methods



# What are the Ways to Sample?

## Traditional sampling: Pros, Cons and Limitations

	Open Sampling Valve	Steam in place valve	Septum sampling	Aseptic Connectors	Tube Welding
					
	<p><b>Large number of samples</b> Low cost per sample</p>	<p>Large number of samples <b>Closed sampling</b> Low cost per sample</p>	<p><b>Large number of samples</b> Low cost per sample</p>	<p>Flexible &amp; Reliable Aseptic sampling <b>Safe and disposable</b></p>	<p>Large number of samples <b>Aseptic sampling</b> Flexible</p>
	<p>Dead-leg <b>Loss of product</b> (flush) Open sampling Impossible to sterilize</p>	<p>Complex operation (SIP) Safety hazard (heat) <b>Risk of sample dilution</b> Container limitations</p>	<p>Not steam sterilizable Safety hazard (needles) <b>Sample volume limitation</b></p>	<p><b>Extra cost</b> Potential dead-leg (tubing) Limited disconnection</p>	<p><b>Waste of product</b> Requires utility Piece of hardware</p>
	<p><b>High risk of contamination</b> Operator &amp; Process safety Sample representativeness</p>	<p>Operator training &amp; safety <b>Sample representativeness</b></p>	<p><b>High risk of contamination</b> Operator safety <b>Sample representativeness</b></p>	<p><b>Operator training</b></p>	<p><b>Requires maintenance</b> Operator training</p>

# Complexity and risks of traditional sampling methods

- FDA Warning letters (483) issued in 2015
- 104 – Procedures designed to prevent microbiological contamination of sterile drug products not established, written or followed
- 24 – Representative samples not obtained

## Microbial contaminations during manufacturing

- Increases risk
- Results in enormous cost
- Requires complicated investigations to prevent reoccurrence



**Cost of sampling VERSUS cost of a shut down**

# Biologics in-process contamination



30  
Percent

*Percent of process deviations  
caused by contamination\**

*Length of time to complete an  
investigation*

1-6  
Months

*Operations cost*

1-14  
Million Euro

\*Source Langer 2013, Wiebe 2014

# Biologics in-process contamination



30

Percent

1-6

Months

1-14

Million Euro

## *Impact*


- ◆ Productivity losses
- ◆ Material replacement costs
- ◆ Batch loss
- ◆ Interruption of product supply
- ◆ Delay in clinical development

# Key considerations for Closed & Disposable sampling options

# Considerations for adopting single-use Aseptic sampling

Situations	Implications
Contamination of samples (out of in process specifications)	<ul style="list-style-type: none"> <li>• Product batch on-hold</li> <li>• Internal investigations</li> <li>• Additional tests on final product</li> <li>• Conflict between manufacturing and QAQC</li> <li>• Risk of remarks from regulatory inspectors</li> </ul>
Contamination of process	
Complex sampling procedure	<ul style="list-style-type: none"> <li>• Risk of false results</li> <li>• Frequent operator training</li> <li>• Unhappy operators</li> <li>• Internal investigations</li> </ul>

# Considerations for adopting single-use Aseptic sampling

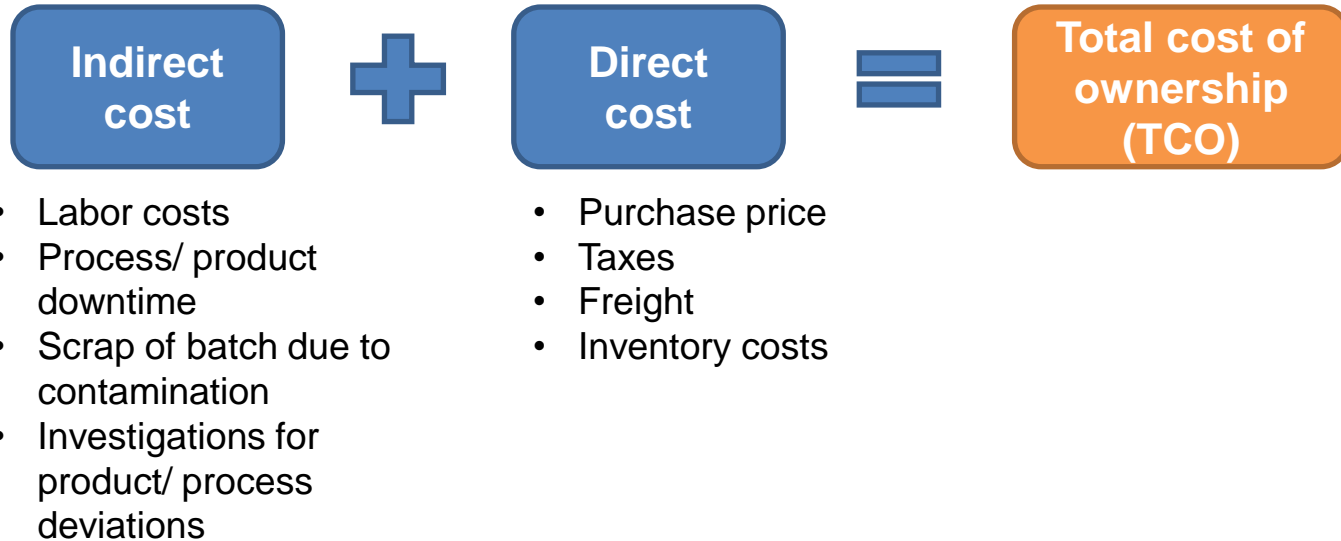

# Self-assessment: Am I suitable to adopt single-use Aseptic sampling?

Considerations	Strong NO → Strong YES			Considerations	Strong NO → Strong YES		
	1	2	3		1	2	3
It is difficult to implement changes to my infrastructure (e.g. piping).				There is high risk of bio-contamination in my process.			
There are challenges to conduct cleaning validation.				My current sampling method has risk of cross-contamination.			
My operators are very familiar with plastic components in the process.				I need to collect samples anywhere.			
I have limited process time to prepare sampling assemblies (traditional).				I need to collect large numbers of samples.			
I have limited cycle time. I need quick turnaround between batches.				I require a varied range of sampling containers.			
I am concerned about operator safety.				My process is sensitive to price.			
I have limited resources for operator training.							



# Sterile sampling value assessment

## Case study example

### Background

Replace standard open sampling with NovaSeptum® sterile sampling assembly

### Outcome



# 10%

## Overall cost reduction

### Key facts

# 62%

Reduced labor hours

# 35%

Reduced product loss during sampling

# 80%

Reduced deviation costs

# 400

Autoclave cycles eliminated

[NovaSeptum Total Cost of Ownership](#)  
[Youtube Video](#)

# Conclusions

- There are complexity and heightened risks in using traditional sampling methods.
- Regulatory recommendations are in place and single-use aseptic sampling could help meet/exceed those corresponding needs.
- There are several key considerations to design/optimize your sampling plan.

# Acknowledgements

Mark Antoine Kaag, Global Product Manager, France

Janmeet Anant, Global Regulatory Advocate, USA

# Thank You

[Somasundaram.g@merckgroup.com](mailto:Somasundaram.g@merckgroup.com)

